

# Analysis Software Installation

## FROM WEBSITE/EMAIL

- **Save** the **QFT\_vX.XX.X\_setup.zip** file (where **X.XX.X** denotes the version of software being installed) to an appropriate location on the computer's hard drive.
- Using **My Computer**, locate the **QFT\_vX.XX.X\_setup.zip** file and double-click on it. This will open the program Winzip®, which can be used to unzip the QFT Analysis Software **installation files**.
- Unzip the installation files to an appropriate location on the computer's hard drive—for example, the **temp** folder.
- Using My Computer, locate and run the file **QuantIFERON\_Startup.exe**. An installation screen, as outlined below, will appear.

## FROM SELF-INSTALLING CD-ROM

- Insert the QuantIFERON-TB Gold In-Tube Analysis Software CD-ROM into the CD-ROM drive. An installation screen will be automatically displayed.
- Select the option **Run QFT vX.XX.X Installation**. Follow the prompts to install the software and support files onto the computer. A folder called **Program Files\QuantIFERON** is created for this purpose.  
Shortcuts to the Analysis Software are created on the desktop and in the Start Menu.

Alternatively, from the **Start Menu** select **Run** and then **Browse** to locate the Analysis Software .zip file. Ensure that **All Files** is chosen in the **Files of Type** field, then select **OK**.

If not currently installed, the program Winzip can be obtained from the website [www.winzip.com](http://www.winzip.com)

The software and support files can also be accessed directly from the **CD-ROM** by selecting the other options on the installation screen.

For more information, please contact the Cellestis office nearest you or visit [www.cellestis.com](http://www.cellestis.com).

**North America / South America**  
Cellestis Inc.  
Email: [customer.service@cellestis.com](mailto:customer.service@cellestis.com)  
Fax: +1 661 775 7479  
Tel: +1 661 775 7480 (outside USA)  
Toll free: 800 519 4627 (USA only)

**Europe / Middle East / Africa**  
Cellestis GmbH  
Email: [europe@cellestis.com](mailto:europe@cellestis.com)  
Fax: +49 6151 428 59 110  
Tel: +49 6151 428 59 0

**Asia / Oceania**  
Cellestis International  
Email: [quantiferon@cellestis.com](mailto:quantiferon@cellestis.com)  
Fax: +61 3 9571 3544  
Tel: +61 3 9571 3500